

DISCOVERY AND ENGINEERING OF ANTI-CD3 HEAVY CHAIN-ONLY ANTIBODIES FOR USE IN T-CELL ENGAGING THERAPEUTICS

ADIMAB

Paul Khalifé*, Hannah Watkins*, Elizabeth McGurk, Cameron Henkel, Todd Boland, Hailey Heston, Whitney Renaud, Christin Strong, Kaleigh Canfield, John Martel, Chanita Sandberg, Rebecca Niles, Liz Parker, Jingfu Zhao, Carissa Metzger, Xiaojun Lu, Irina Burnina, Asparouh Lilov, Allie Karakosta, Hannah Green, Kelsie Leary, Melanie Sinclair, Shu Lin, Mike Brown, Jessica Dawson, Cory Ahonen, Andrew Avery, Mike Battles, Dan Walsh, Garrett Rappazzo, Nicole Amagliani, Lauryn Taylor, Sarat Pudi, Robert Pejchal, Beth Sharkey, Morgan Morrill, and Noel Pauli

INTRODUCTION TO T CELL ENGAGERS

T cell engagers (TCEs) utilizing CD3 are an increasingly validated class of multispecific antibodies that have shown promise in treating oncologic conditions. Through pairing an arm targeting CD3 on T cells with an arm targeting a tumor associated antigen, the TCE forces the T cell and tumor cell into close proximity, triggering T cell activation leading to tumor cell killing¹.

- Eight TCEs are currently FDA approved- six since 2022¹
- TCEs can be engineered in multiple bi- and multispecific formats. Activation level can be influenced by format¹

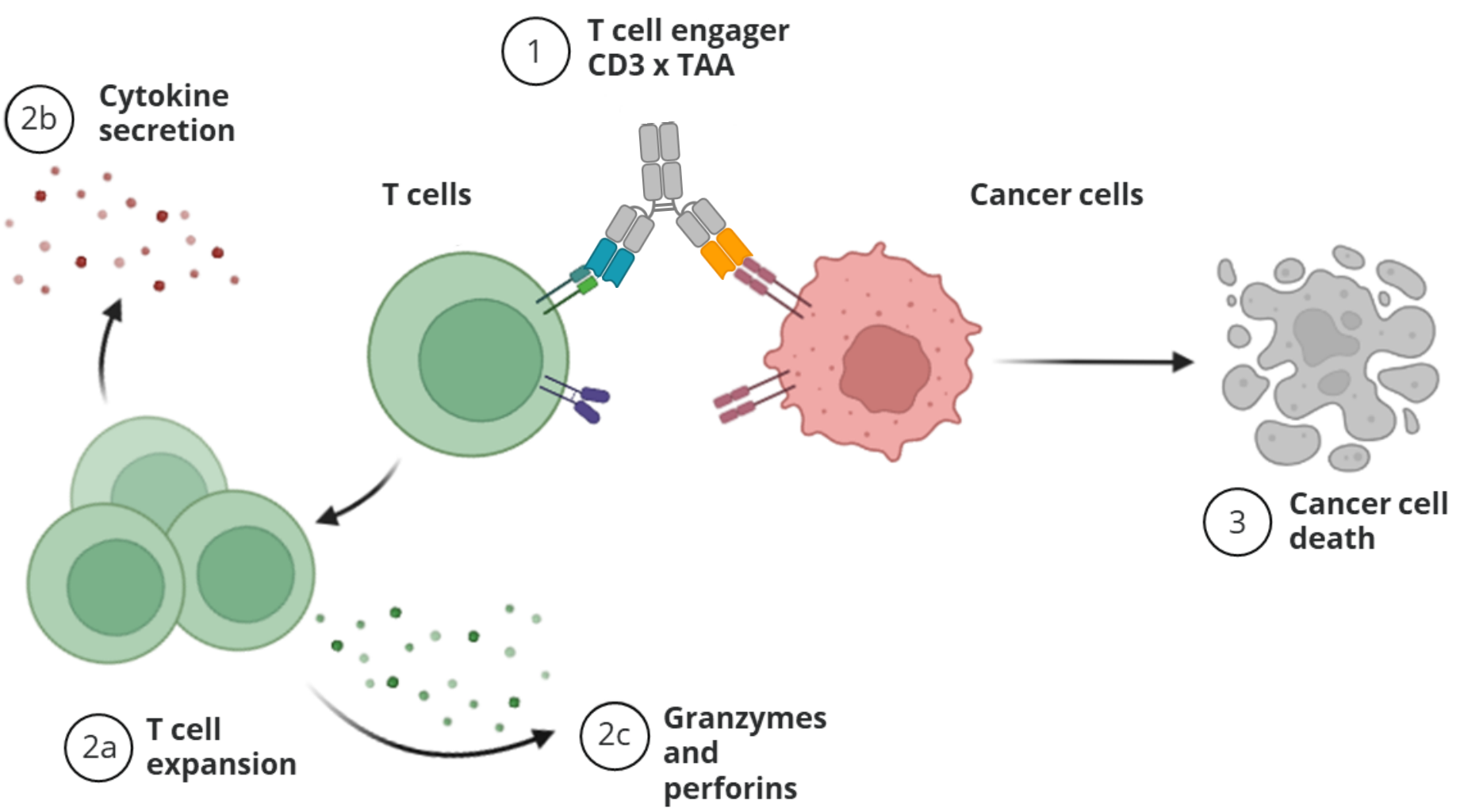


Figure 1. Summary of T cell engager mechanism of action.

ADIMAB α-CD3 CONVENTIONAL IgG LINEAGE: HUMAN x CYNO CROSS-REACTIVITY AND TUNABLE POTENCY

In vivo murine discovery followed by humanization and yeast-based optimization generated a broad affinity α-CD3 panel with excellent developability properties.

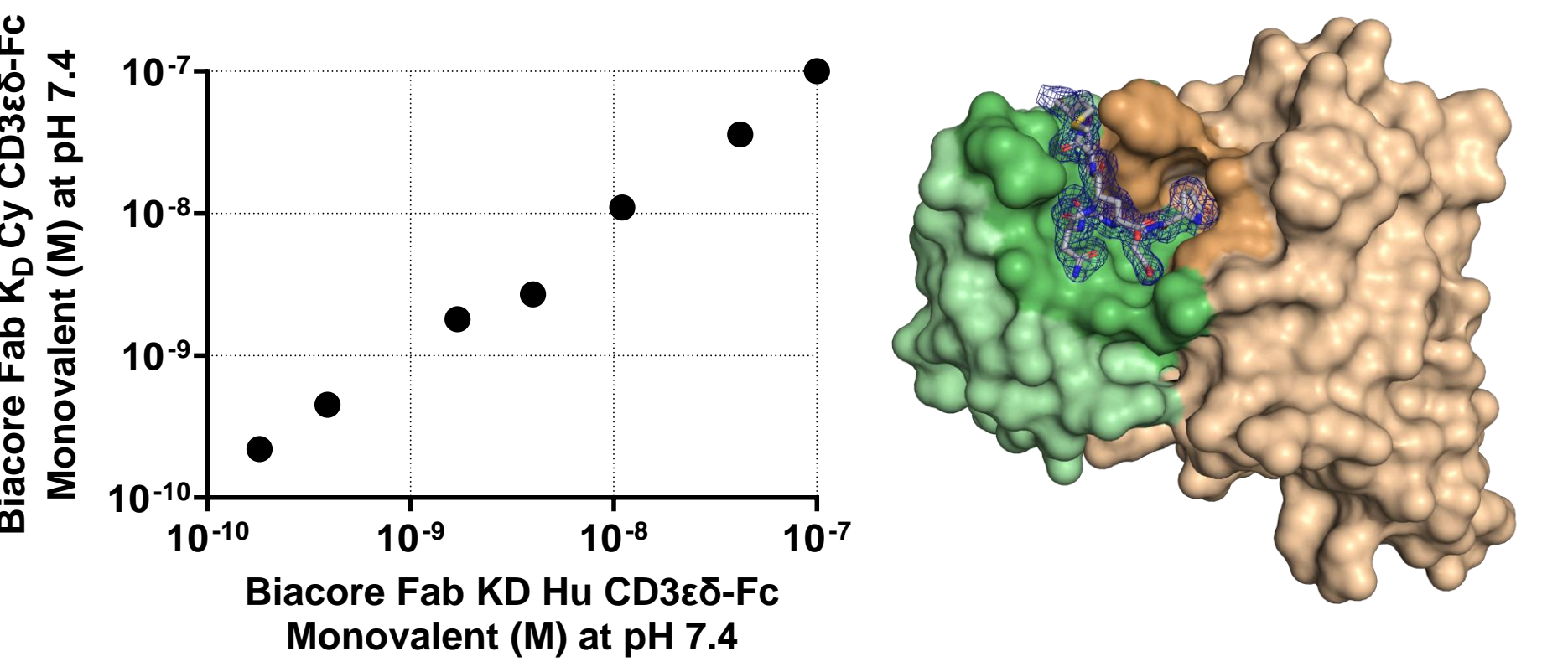


Figure 2. Adimab's α-CD3 antibody lineage has human-cyno cross-reactivity across a broad range of affinities and binds to the N-terminal portion of CD3ε².

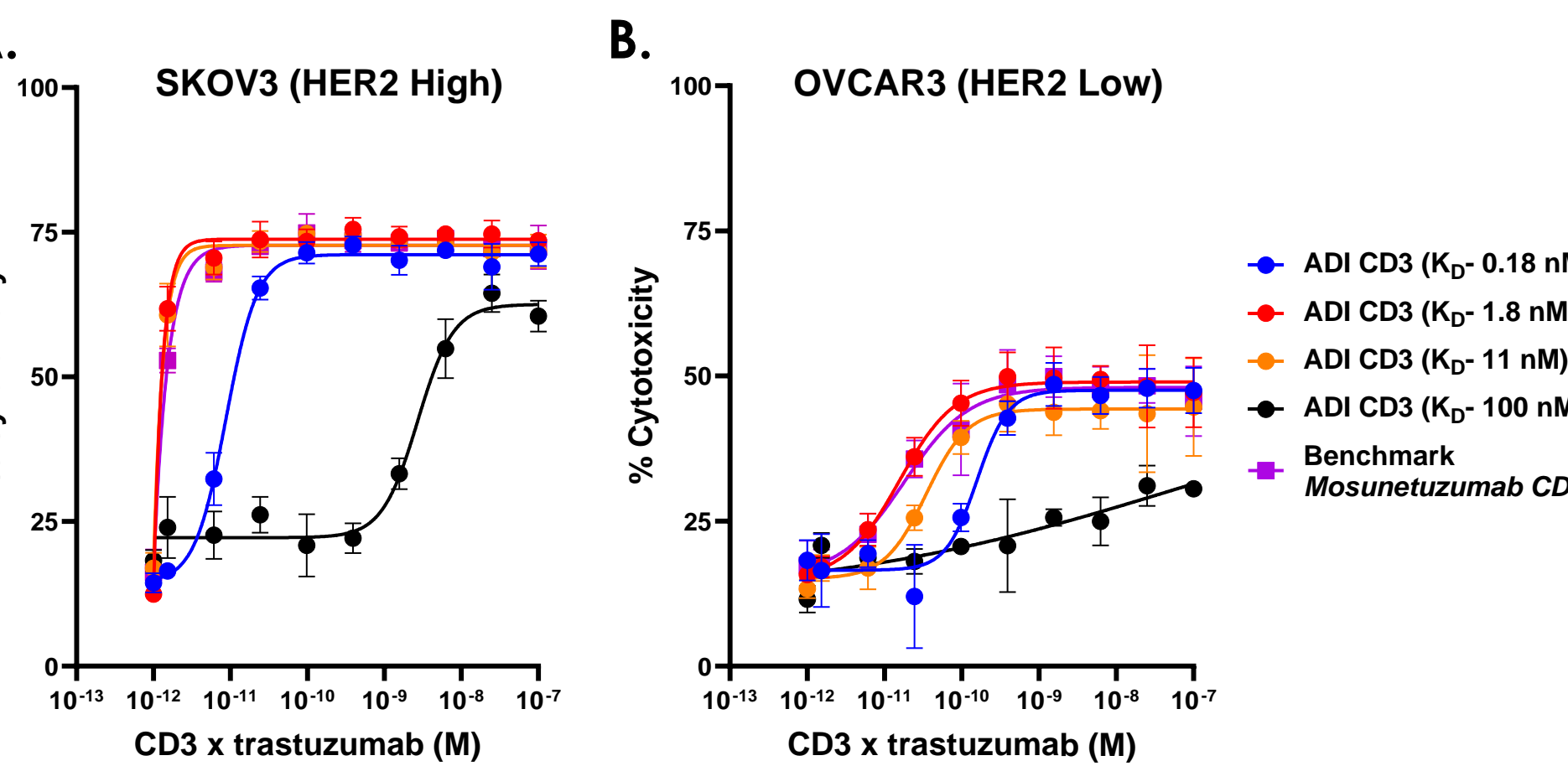


Figure 3. CD3 x HER2 bsAbs support potent Redirected T Cell Cytotoxicity (RTCC) on A. SKOV3 (HER2 high) and B. OVCAR3 (HER2 low) cell lines. The monovalent K_D for CD3 for each bispecific IgG is indicated in the legend. α-HER2 Fv is trastuzumab. Effector:target (E:T) 5:1, 24-hour time point shown.

EXPANDING ADIMAB'S CD3 BINDERS: HCAB DISCOVERY

Camelids naturally produce two classes of antibody: conventional (HC+LC) and HCABs (HC-only)³. Absence of a light chain can simplify bi- or multispecific development⁴. Adimab's yeast-based immune library platform facilitates discovery of llama-derived HCABs.

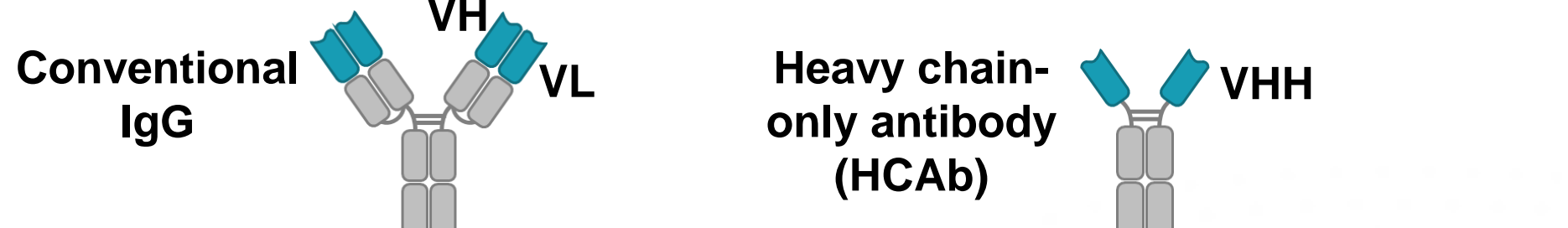


Figure 4. Comparison of conventional IgG to HCAB

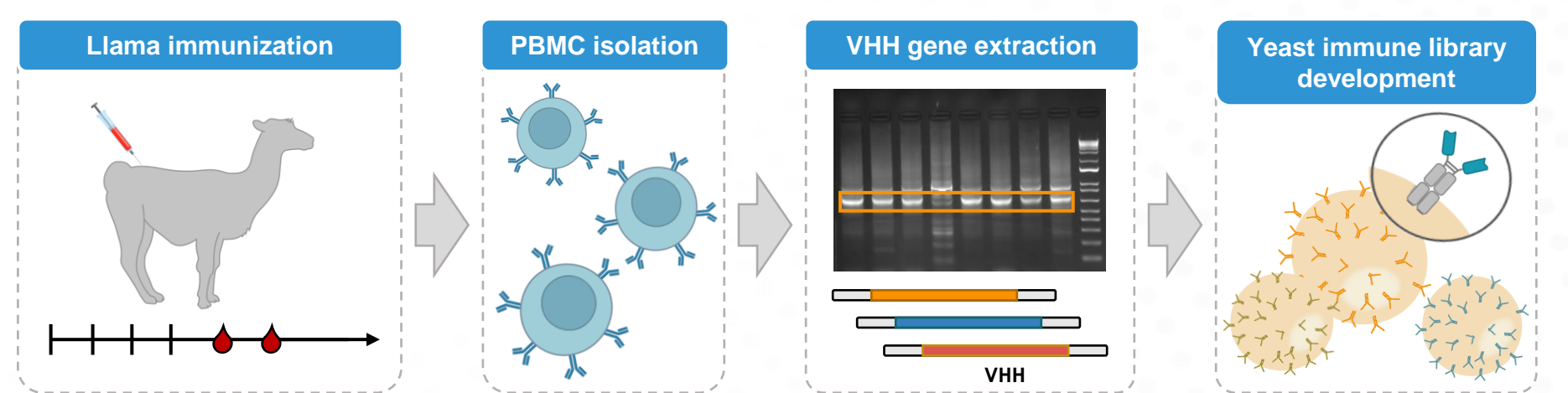


Figure 5. Building of α-CD3 immune libraries using Adimab platform

IMMUNE LIBRARY SELECTIONS DISCOVERED HUNDREDS OF α-CD3 HCABS

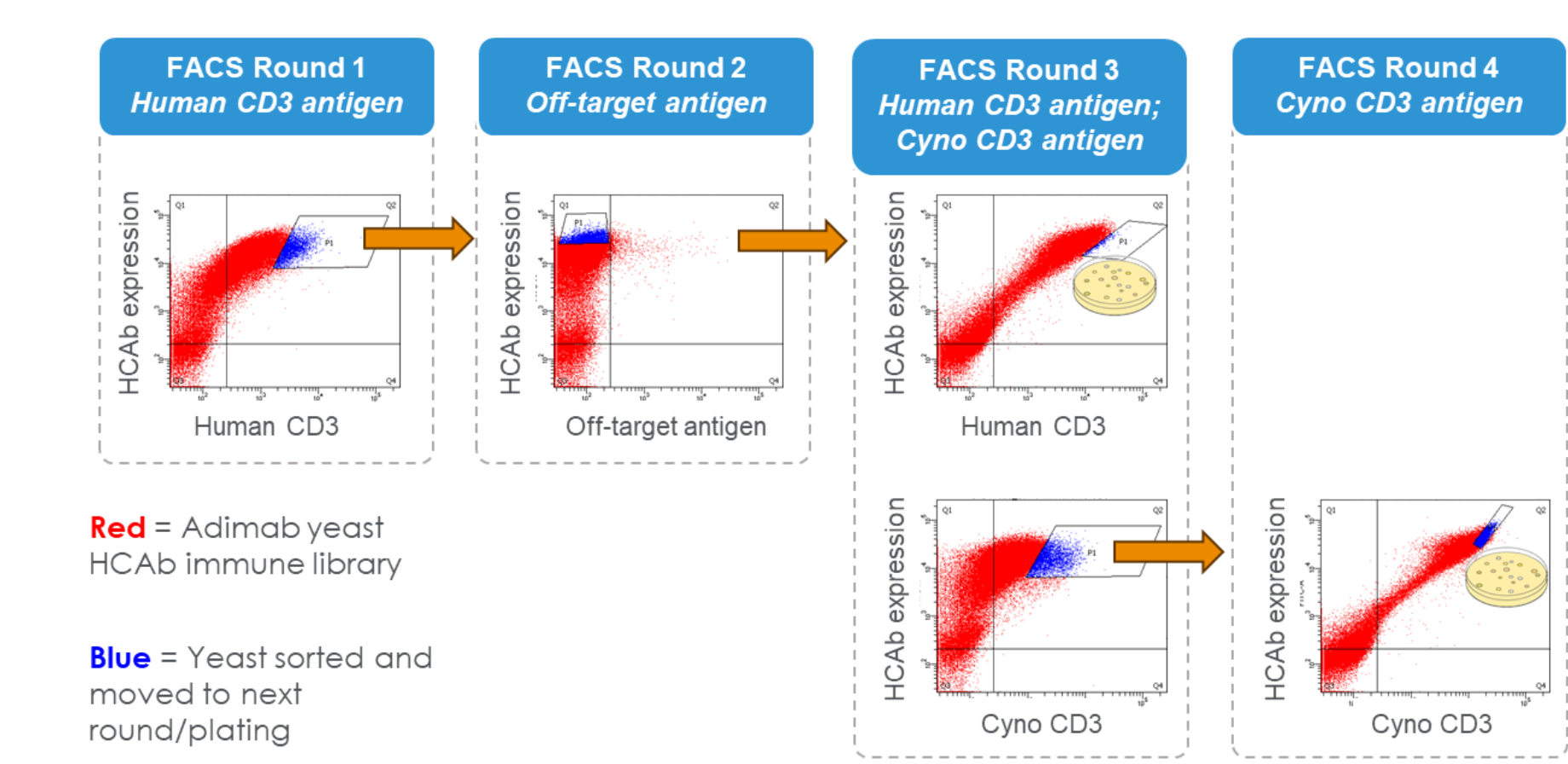


Figure 6. Multiple rounds of selections were performed on the Adimab yeast immune libraries to isolate α-CD3 HCABs with desirable properties. Selections were performed under conditions to push for cyno cross-reactivity, wide affinity range, and favorable developability. Note: figure displays only one representative selection.

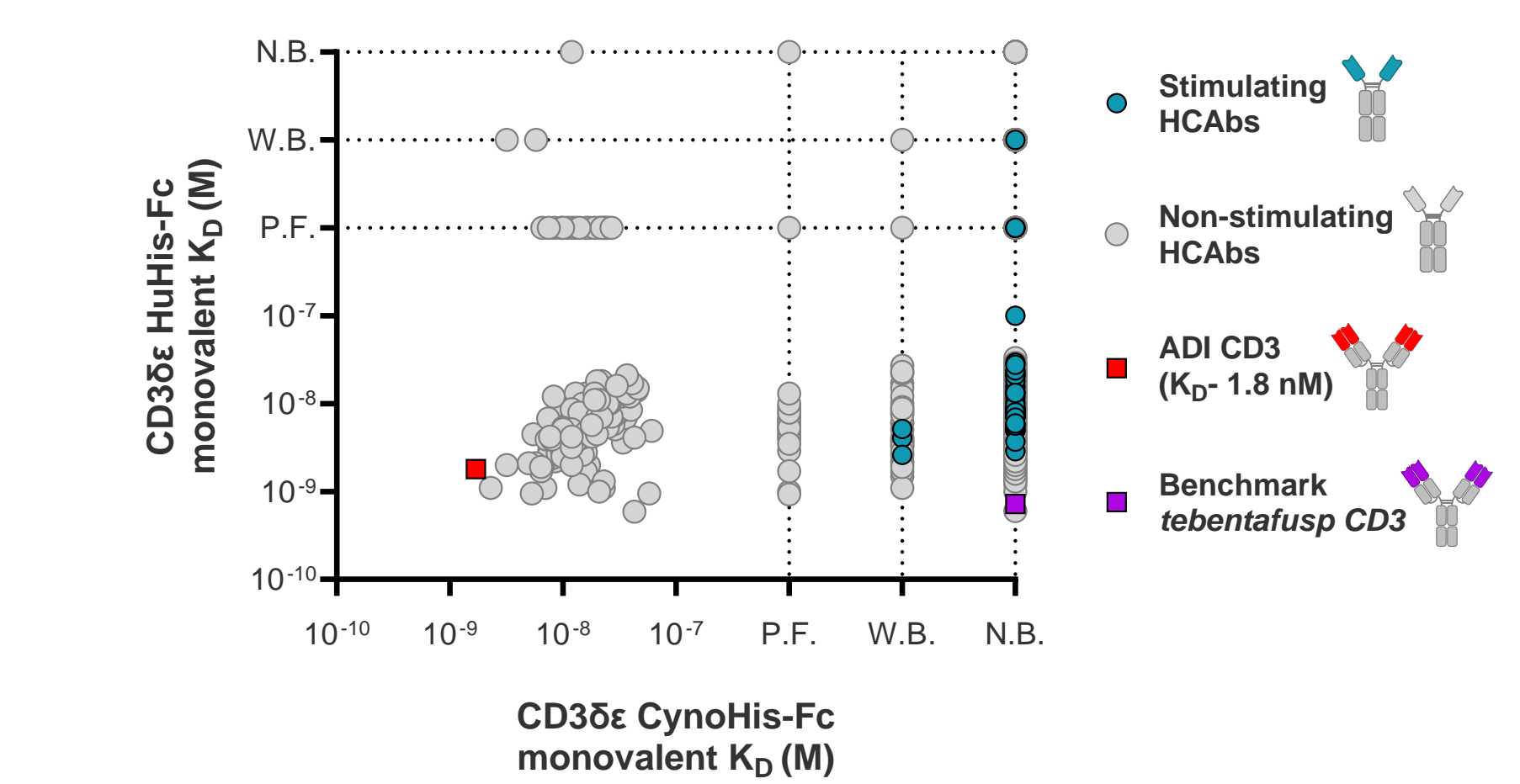


Figure 7. Isolated HCABs screened for affinity and activity in CD69 upregulation T cell stimulation assay. While many human x cyno α-CD3 binding HCABs were isolated, stimulating HCABs exhibited only human-specific binding. P.F. = poor fit; W.B. = weak binding; N.B. = no binding.

α-CD3 T-CELL ENGAGERS SHOW ACTIVITY IN FUNCTIONAL ASSAYS AND DISPLAY FAVORABLE DEVELOPABILITY

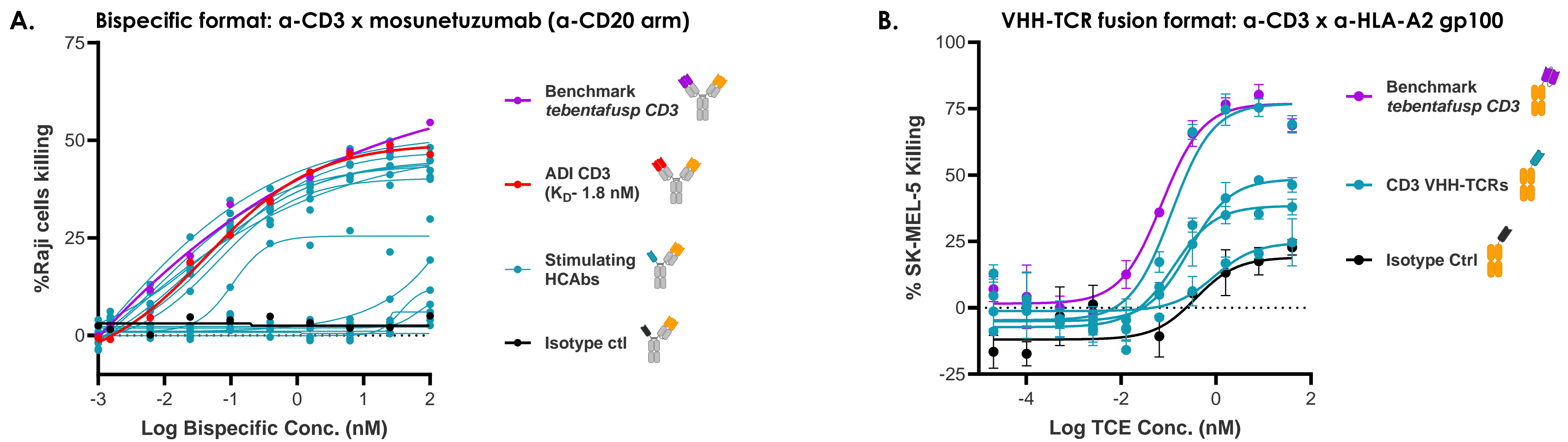


Figure 8. Stimulating HCABs were downselected, reformatted in Chain Exchange vectors, and produced in two different TCE formats: as α-CD3 HCAB x α-CD20 bispecific antibodies (bsAbs) and as VHH-T-cell receptor fusions (VHH-TCRs). α-CD3 x α-CD20 bsAbs evaluated in RTCC assays resulted in cytotoxicity against target Raji (CD20+) cells (A.). VHH-TCRs evaluated in RTCC assays resulted in cytotoxicity against melanoma (SK-MEL-5) cells (B.).

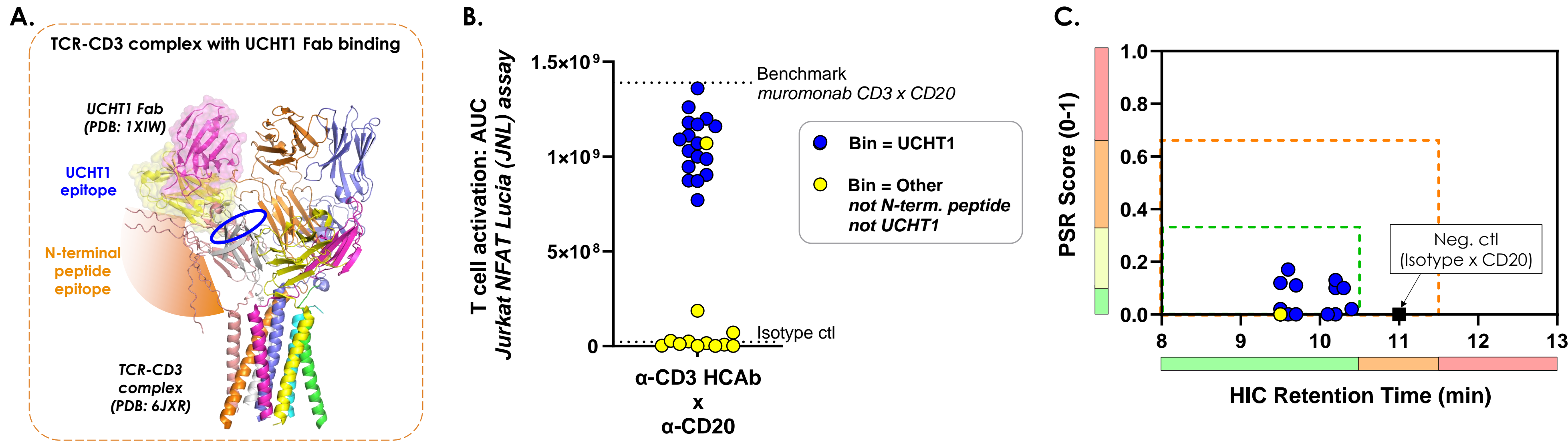


Figure 9. A. TCR-CD3 complex shown with an overlay of UCHT1 (basis for the CD3 arm of tebentafusp); schematic of N-terminal peptide also shown (target of the CD3 arm of mosunetuzumab). B. The majority of Adimab-discovered functional HCABs competed with UCHT1, and none bound the N-terminal peptide. C. All clones with functional activity in the JNL and RTCC assays displayed favorable developability with low PSR and HIC retention times⁵.

HUMANIZED α-CD3 HCABS MAINTAIN ABILITY TO STIMULATE T CELLS

Six functional HCABs, spanning six lineages and two bins, were selected for humanization. A panel of humanized constructs were designed for each parent, retaining the llama CDRs and select amino acids within the frameworks.

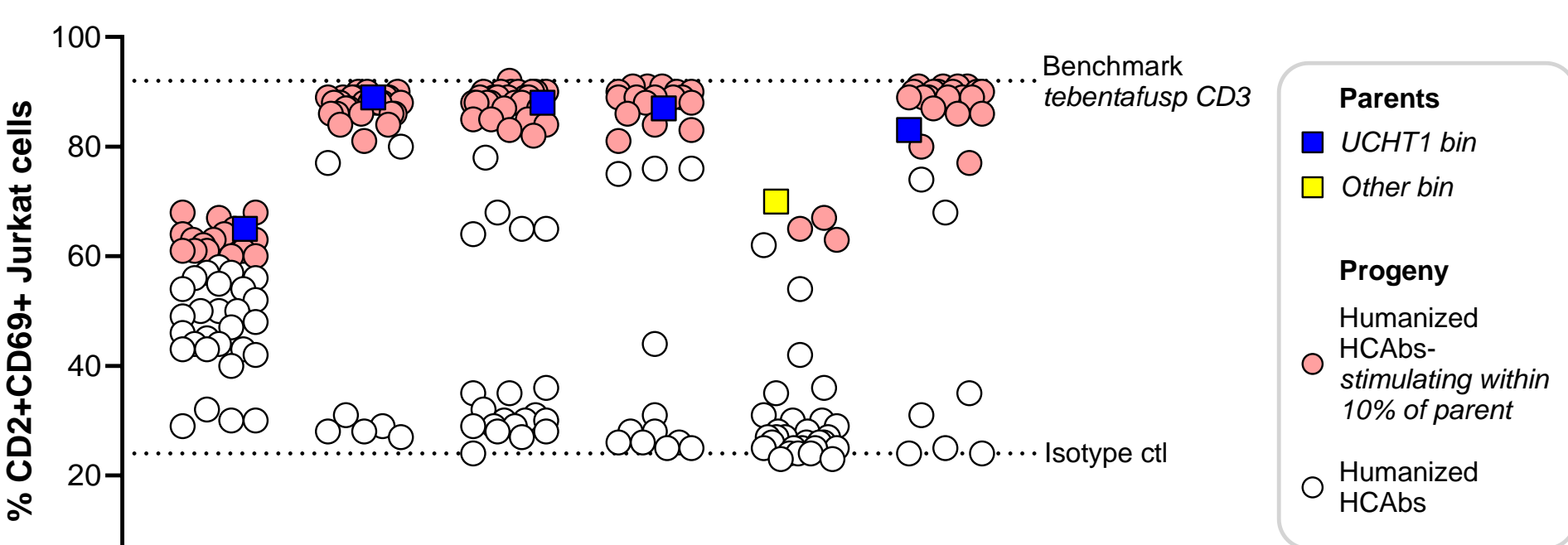


Figure 10. Humanized HCABs were screened in a CD69 upregulation T cell stimulation assay. All parents had progeny that stimulated within 10% of parent.

ACKNOWLEDGEMENTS

Thanks to the following teams and cores at Adimab for their contributions to this work: in vivo discovery, yeast and mammalian protein production, protein analytics, sequencing and molecular, computational biology, product engineering, legal, alliance management, and business development.

SUMMARY

Heavy chain-only antibodies (HCABs) offer promise in reducing the complexity of multispecific therapeutics. Here, we describe the discovery and engineering of a panel of high-affinity, CD3-specific HCABs from immunized llamas, and their validation for use in a T cell engaging bispecific format. Further validation and optimization are in progress. The resulting set of HCABs, and future improvements to the panel, can be accessed as part of Adimab's non-exclusive TCE offering.

REFERENCES

- Center for Drug Evaluation and Research. Bispecific Antibodies: An Area of Research and Clinical Applications, FDA, 14 Feb. 2024, www.fda.gov/drugs/news-events-human-drugs/bispecific-antibodies-area-research-and-clinical-applications.
- Liu, CY, et al. 2023. Structure-based engineering of a novel CD3ε-targeting antibody for reduced polyreactivity. MABS 15(1): 2189974. PMID: 36991534.
- Hamers-Casterman, C. T. S. G., et al. (1993). Naturally occurring antibodies devoid of light chains. Nature, 363(6428), 446-448. PMID: 8502296.
- Wesolowski, J. et al. (2009). Single domain antibodies: promising experimental and therapeutic tools in infection and immunity. Medical microbiology and immunology, 198, 157-174. PMID: 19529959
- Jain, et al. (2017). Biophysical properties of the clinical-stage antibody landscape. Proceedings of the National Academy of Sciences, 114(5), 944-949. PMID: 28096333