Pairing in vivo diversities with a yeast-based platform for integral membrane protein-specific antibody discovery



Noel Pauli, Jiannan Li, Hannah Watkins, Kaleigh Canfield, Todd Boland, Cory Ahonen, Rebecca Niles, Robert Pejchal, and Eric Krauland (Adimab LLC, Lebanon, NH, USA)

BACKGROUND

Integral membrane proteins continue to be a hurdle for antibody therapeutic discovery. By combining a yeast-based, B-cell platform with diversities derived from humanized transgenic murine or wildtype mice, and llamas, we have developed a high-throughput methodology for the discovery of membrane-obligate-target (MOT) specific antibodies. In the absence of soluble recombinant antigen, this platform enables the discovery of large panels of clonally-diverse, high-affinity, target-specific IgGs or Heavy Chain-only Antibodies (HCAbs).

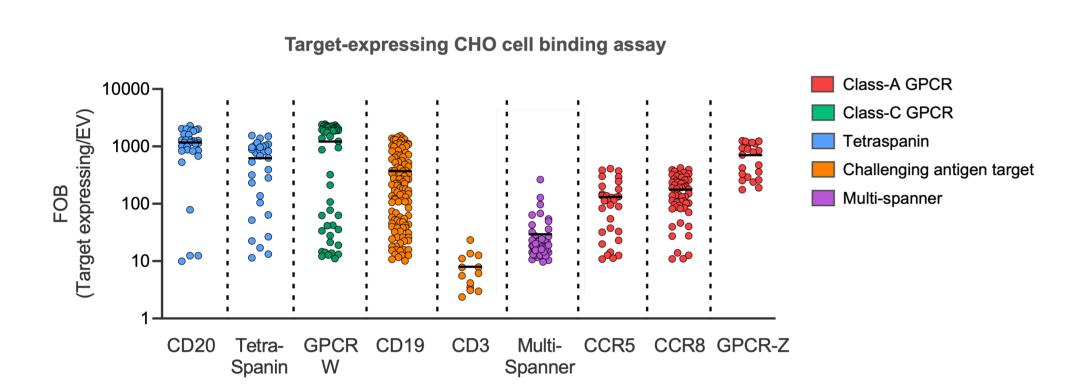


Figure 1: Discovery of mAbs against different types of MOTs, including GPCRs and tetraspanins.

Summation of multiple membrane-obligate projects where the Y-axis displays fold-over-background (FOB) binding signal of antibodies used to stain target-expressing CHO cells versus Empty Vector (EV) CHO cells at 100 nM staining concentration

METHODS

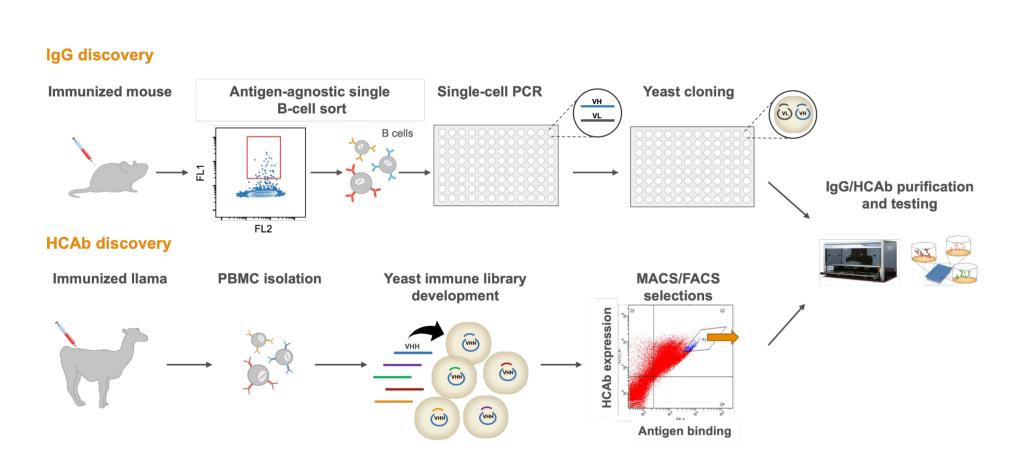


Figure 2: Adimab pairs immunized animal diversities with our engineered yeast-based platform for the rapid discovery of membrane protein-specific mAbs.

RESULTS

CCR8 mAb Discovery and Characterization

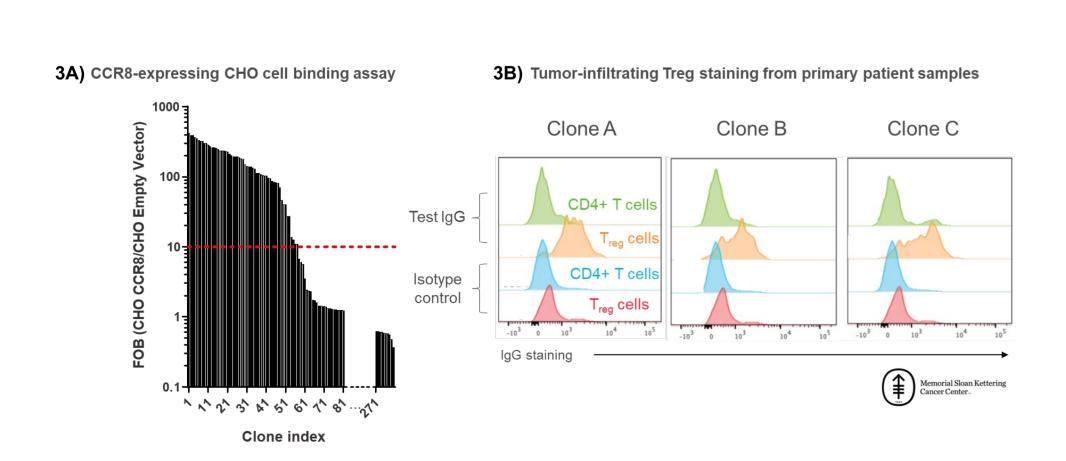


Figure 3: Efficient recovery of many CCR8-specific antibodies using Adimab's MOT platform.

3A) 57 CCR8-specific antibodies recovered following workflow. **3B)** Representative histograms of individual IgG clones stained on tumor infiltrating T cells taken from human donors at 100 nM primary staining.

RESULTS

CCR8 mAb Discovery and Characterization

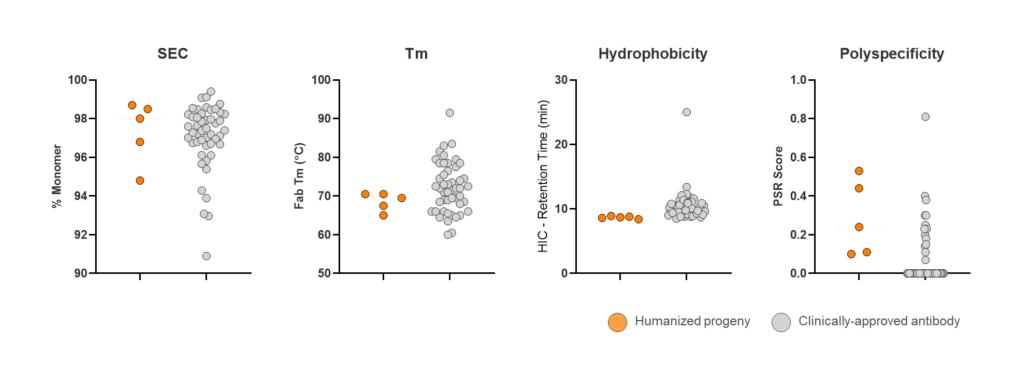


Figure 4: Humanized CCR8-specific IgGs displayed favorable developability profiles.

Results of Size Exclusion Chromatography (SEC), melting temperatures (Tm), Hydrophobic Interaction Chromatography (HIC), and Poly-Specificity Reagent (PSR) binding, respectively. Antibodies were HEK-produced and compared against clinically-approved antibody developability data (Jain, et al., 2017)

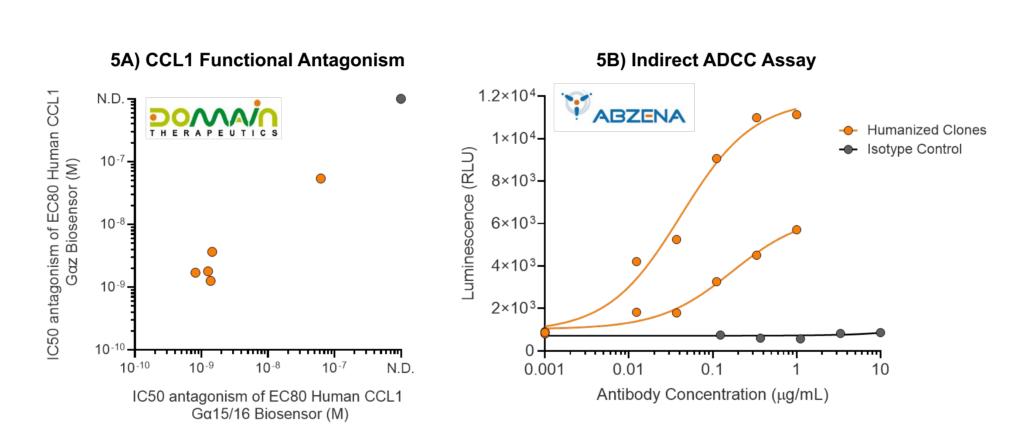


Figure 5: CCR8 mAbs antagonize CCL1 and display ADCC activity.

5A) GPCR Biosensor BRET-based assay which detects CCR8 signaling when in the presence of its natural ligand, CCL1. **5B)** Results of indirect Antibody Dependent Cell-mediated Cytotoxicity (ADCC) assay in which genetically-modified effector cells express luciferase upon antibody-dependent recognition of antigen expressed on target cells

'GPCR-A' Immune Library HCAb Discovery

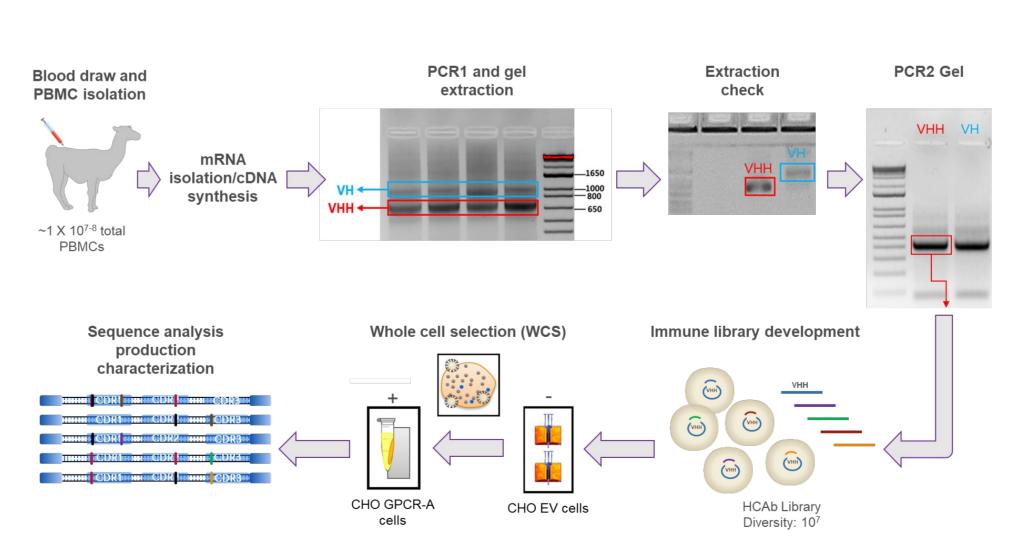


Figure 6: Integration of immunized llama VHH B cell diversities for the generation of yeast immune libraries for selection against MOT "GPCR-A".

Discovered VHH regions are cloned into yeastspecific acceptor vectors that contain a modified hinge fused with the human IgG1 Fc domain to create the HCAb format.

'GPCR-A' Immune Library HCAb Discovery

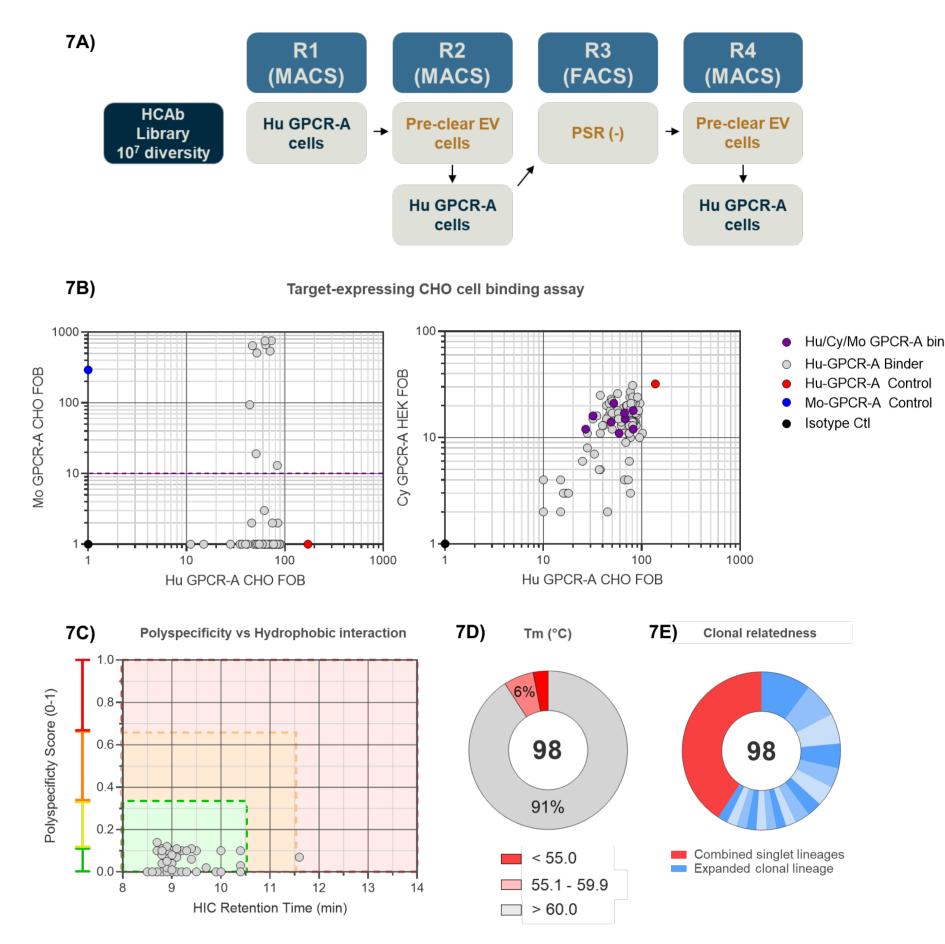


Figure 7: Triple cross-reactive GPCR-A-specific HCAbs were successfully isolated using Adimab's MOT platform.

7A) Four-round immune library selection scheme. **7B**) 98 unique HCAb isolated against GPCR-A were assayed at 100 nM for binding against cell lines expressing Hu, Cy, and Mo GPCR-A. **7C and 7D**) Selection output developability analysis of polyspecificity, HIC retention time, and Tm. **7E**) Clonal lineage analysis comparing sequence relatedness in GPCR-A specific HCAb output

CONCLUSIONS

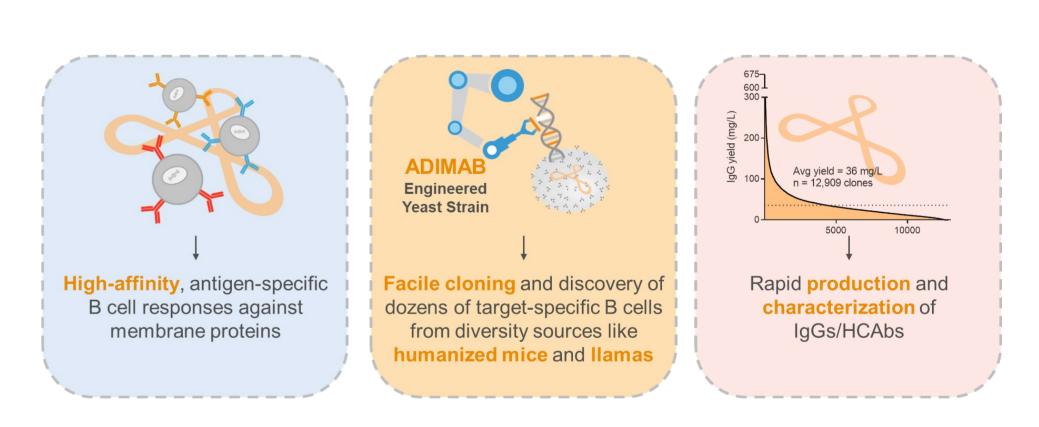


Figure 8: Adimab's yeast-based platform, paired with immunized in vivo diversities, can solve many of the challenges of membrane-protein antibody discovery

REFERENCES

(1) Jain, Tushar et al. "Biophysical properties of the clinical-stage antibody landscape." *PNAS*, vol. 114,5 (2017): 944-949.

ACKNOWLEDGEMENTS

- Memorial Sloan Kettering Cancer Center: George Plitas and Alexander Rudensky
- Domain Therapeutics
- Abzena: Biologics Research Partner